

Sütte Kantitatif Aflatoksin M1

Tespiti için Test kitleri

Aflasensor süt numunelerinde Aflatoksin M1 moleküllerinin kantitatif tespiti için geliştirilmiş hızlı test kitidir.



ÖZELLİKLER

**KANTİTATİF****HIZLI**

10 DAKİKADA SONUÇ.

**HASSAS**

STANDARTLARA UYUMLU TESPİT (AB-ABD).

**GÜVENİLİR VE DAYANIKLI****KULLANICI DOSTU**

SAHADA VEYA LABORATUARDA KOLAY KULLANIM.

**DÜŞÜK MALİYET AVANTAJI**

SONUÇLAR

Readsensor2 AFM1 kontaminasyonun kantitatif tespitini sağlar.



DUYARLILIK

TOKSİN	AB MAKSİMUM KALINTI LİMİTİ	ABD MAKSİMUM KALINTI LİMİTİ
		KIT041 (KANTİTATİF)

Aflatoksin M1 (ppt)

Tespit Aralığı 30-150 ppt

Tespit Aralığı 200-750 ppt

NASIL KULLANILIR




1 Sütü ve reaktifleri karıştırın



2 Kuyucukları inkübatöre yerleştirin



3 Başlığı kapatın ve inkübasyonu başlatın

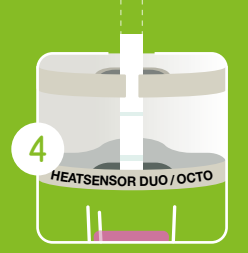

4 Test şeritlerini inkübatöre yerleştirin

3+7 = 10 DAKİKA 40°C DE İNKÜBE EDİN



5 Sonuçları Readsensör2'de okuyun

KANTİTATİF

AflaSensor Quanti

ÜRÜNLER

KIT041v2

AFLASENSOR QUANTI MK 0,05 PPB - 96 PC

KIT078

AFLASENSOR MK 0,5 PPB 96 pc

APP032 HEATSENSOR DUO

APP008 AIRSHIELD DUO

APP052 HEATSENSOR OCTO

APP033 AIRSHIELD OCTO



APP088

READSENSOR 2



unisensor
DIAGNOSTIC ENGINEERING

Liège Science Park
Rue Louis PLescia, 8
B-4102 Ougrée (Liège)
BELGIUM

www.unisensor.be

Phone +32 4 335 30 00
info@unisensor.be

Reseller

mayasan
biotech

Mayasan Gıda San.ve Tic. A.Ş.

Address: Omerli mah. Hadimkoy-Omerli cad. No:18 34555 Arnavutkoy/Istanbul/Turkey

Phone: 0090 212 639 6565 (pbx) • Email: sales@mayasan.com • www.mayasan.com